

Catalog Number: HZ-1151-GMP

## **Data Sheet**



## HumanKine® recombinant human FLT3 Ligand protein- GMP grade

**Animal Component-Free** 

**Human cell expressed** 

Tag-Free


**Endotoxin Free** 

## **Product Description**


Animal-free Recombinant Human FLT3 Ligand is expressed from human 293 cells as a monomeric glycoprotein with an apparent molecular mass of 24 to 30 kDa. This cytokine is produced in a serum-free, chemically defined media. The molecular mass is a narrower range than reported when this cytokine is expressed in insect Sf21 cells (17 to 30 kDa), emphasizing the difference in glycosylation in different expression systems. FLT3 Ligand is a growth factor that regulates proliferation of early hematopoietic cells. FLT3 Ligand binds to cells expressing the tyrosine kinase receptor FLT3. FLT3 Ligand by itself does not stimulate proliferation of early hematopoietic cells, but synergizes with other CSFs and interleukins to induce growth and differentiation.

Alternative Names	FL, Flt3 ligand, Flt3L, FLT3LG, SL cytokine		
Accession Number	P49771		
Source	Human Embryonic Kidney cells (HEK293). HEK293-derived FLT3 Ligand protein		
Species Reactivity	human, mouse		
Adventitious Virus	Adventitious Virus Master Cell Bank tested Negative for Adventitious Viruses		

Specifications					
Test	Method	Specification			
Activity	Dose-dependent proliferation of the human acute myeloid leukemia cell line (OCI-AML5).	0.4-3.0 ng/mL			
Molecular Mass	SDS-PAGE	27 to 34 kDa reduced, 23 to 30 kDa non-reduced, monomer, glycosylated			
Purity	SDS-PAGE	>95%			
Endotoxin	LAL	<0.1 EU/μg			
Mycoplasma	PCR	Not Detected			

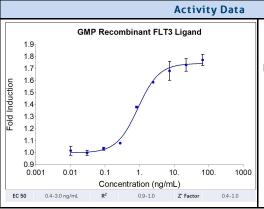

GMP-grade recombinant human FLT3 Ligand (HZ-1151-GMP) stimulates dose-dependent proliferation of the human acute myeloid leukemia cell line (OCI-AML5). Viable cell number was quantitiatively assessed by PrestoBlue Cell Viability Reagent. OCI-AML5 cells were treated with increasing concentrations of recombinant human FLT3 Ligand for 72 hours. The EC50 was determined using a 4- parameter non-linear regression model. Activity

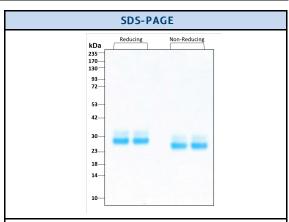

Purity of GMP-grade recombinant human FLT3 Ligand was determined by SDS- polyacrylamide gel electrophoresis. The protein was resolved in an SDS- polyacrylamide gel in reducing and non-reducing conditions and stained using Coomassie blue.

www.ptglab.com

Document #: FR-QA118-101 Rev 0
Data Sheet Version #: 1

Proteintech Group, Inc.

5500 Pearl Street, Suite 400 Rosemont, IL 60612 t: 1-888-478-4522; f: 1-312-455-8408 Email: proteintech@ptglab.com

Preparation				
Shipping Temperature				
Formulation	n 1x PBS, See Certificate of Analysis for details			
Reconstitution Briefly centrifuge the vial before opening. It is recommended to reconstitute the protein to 0.2 mg/mL in sterile water conta endotoxin-free recombinant human serum albumin (HSA). Gently swirl or tap vial to mix.				

	Product Form	Temperature Conditions	Storage Time (From Date of Receipt)
	Lyophilized	-20°C to -80°C	Until Expiry Date
Stability and Storage	Lyophilized	Room Temperature	2 weeks
-	Reconstituted as per CofA	-20°C to -80°C	6 months
	Reconstituted as per CofA	4°C	1 week
	Avoid repeated freeze-thaw cycles.		

## Proteintech GMP Quality Policy HumanKine® GMP Proteins

Invitro recombinant protein production can be prone to variability due to various reasons ranging from quality of raw materials to inconsistency in the process. Therefore, HumanKine®, a proteintech brand's GMP proteins are produced and tested under an ISO 13485 certified quality management system in a clean room facility. Proteintech manufactures the GMP HumanKine® products according to the applicable sections in the following documents: USP Chapter 1043 (Ancillary Materials for Cell, Gene, and Tissue-Engineered Products, USP Chapter 92 (Growth Factors and Cytokines Used in Cell Therapy Manufacturing), WHO TRS, No. 822, 1992 Annex 1 (Good Manufacturing Practices for Biological Products), Ph. Eur. General Chapter 5.2.12, and EudraLex – Volume 4 – Part IV (Guidelines on GMP specific to ATMPs). Proteintech strives to achieve the utmost quality GMP raw material ensuring all applicable guidelines are taken into consideration.

The QMS is built to provide our customers with consistent and pure product delivered by documented processes, qualified personnel, validated processes, qualified equipment, qualified vendors, and a stringent final product release process. Although the final product release process is important, Proteintech performs in-process QC steps after each major manufacturing stage. Production records and facilities may be available for an inspection by approved personnel.

Our GMP policy covers all the aspects of production; from raw materials, facility, equipment, and Instruments to training and personal hygiene of staff. It also guarantees that the process is explicit, validated and well documented for transparency and traceability.

www.ptglab.com

Document #: FR-QA118-101 Rev 0 Data Sheet Version #: 1 Proteintech Group, Inc.

5500 Pearl Street, Suite 400 Rosemont, IL 60612 t: 1-888-478-4522; f: 1-312-455-8408 Email: proteintech@ptglab.com