

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

Introduction

The ChromoTek Nano-CaptureLigand™ human IgG/rabbit IgG, Fc-specific VHH, biotinylated is used for the site-directed and specific immobilization of the Fc-fragment of human IgG and rabbit IgG in biosensor and ELISA assays. It captures non-biotinylated human IgG or rabbit IgG antibodies or Fc-fragments to streptavidin/avidin.

Nano-CaptureLigand human IgG/rabbit IgG, Fc-specific VHH, biotinylated comprises a monoclonal biotinylated VHH/ Nanobody. The product belongs to the Nano-CaptureLigands™ family.

Properties

Description	Monovalent, recombinant single domain antibody for the immobilization of human IgG and rabbit IgG: alpaca monoclonal Nanobody, Fc-specific, biotinylated
Product Type	Capture Nanobody (VHH)
Applications	Immobilization of human and rabbit antibodies on avidin and streptavidin surfaces for Bio-Layer Interferometry (BLI), Surface Plasmon Resonance (SPR) and ELISA
Target / Specificity	Fc-fragment of human IgG1, IgG2, IgG3, and IgG4 and rabbit IgG
Cross-reactivity	No cross-reactivity to human IgA, IgE, IgM; mouse IgG1, IgG2a, Ig2b, Ig2c, IgG3; rat IgG1, IgG2a, IgG2b, IgG2c; sera from goat, guinea pig, rat, sheep Cross-reactivity to macaque (cynomolgus monkey) serum
Affinity (Kd) of monovalent (1:1) binding mode	Human IgG1: 1.4 nM; Human IgG2: 0.4 nM; Human IgG3: 9.0 nM; Human IgG4: 0.2 nM; Rabbit IgG: 0.2 nM Apparent affinity may be higher for full IgGs due to avidity effects (1 antibody captured by 2 Nanobodies).
Concentration	1 g/L (71 µM)
Conjugate	Biotin
Degree of biotinylation	On average 1-2 biotin molecules per Nanobody
Format	Alpaca single domain antibody, monovalent
Host	Alpaca-derived, recombinantly produced in bacteria
Clonality	Monoclonal

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

Clone	CTK0101 (VHH0238)
RRID	AB_2848185
Synonyms	Alpaca single domain antibody, VHH, Nanobody, binding domain of single domain antibody, Nano-antibody
Validation	Application validated for ELISA and BLI (FortéBio Octet® systems) Determination of cross-reactivity, subclass specificity, sequence, affinity, and melting temperature
Purity	Recombinantly expressed and purified via His-tag
Form	Buffered aqueous solution
Storage buffer	25 mM TAPS pH 8.5, 500 mM NaCl, 5 mM EDTA, Preservative: 0.09 % sodium azide
Storage conditions	Upon receipt store at +4°C/+40°F. <i>Optional:</i> Aliquot upon arrival and store at -20°C/-4°F
Stability	Stable for 1 year at +4°C/+40°F
Shipment	Shipped at ambient temperature

Product sizes

Product	Product code	Size
Nano-CaptureLigand™ human IgG/rabbit IgG, Fc-specific VHH, biotinylated	shurbGB-1-10	10 µL
	shurbGB-1-100	100 µL

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

Suggested buffer compositions

Recommended buffers for BLI

Buffer	Composition
1x Kinetics buffer	PBS, 0.01 % (m/v) BSA, 0.002% (v/v) Tween-20
Regeneration buffer	0.01 M glycine, pH 2

Recommended buffers for ELISA

Buffer	Composition
Blocking buffer	PBS, 3 % BSA
Dilution buffer	PBS, 0.05 % Tween-20, 0.5 % BSA
Wash buffer for avidin coating	PBS
Wash buffer	PBS, 0.05 % Tween-20

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

Bio-Layer Interferometry (BLI) protocol

General notes

- The following protocol was designed for the use with a FortéBio Octet Red96e system. Other BLI systems or your specific research question may require optimization of particular parameters.
- Use the recommended materials or their equivalents.
- Set up all samples in a black 96-well microplate (e.g. Greiner Microplate 96 well, PP, flat-bottom, black, #655209) at room temperature. Use 200 µL per well.
- Nano-CaptureLigands are highly compatible with avidin or streptavidin sensors (e.g. FortéBio Streptavidin (SA) Biosensors, #18-5019) and FortéBio Octet® and BLItz® systems.
- Run all experiments at +30°C, a shaking speed of 1000 rpm and a recording rate of 5 Hz.
- Dilute all samples in 1x Kinetics buffer. *Optional:* Use 10x Kinetics buffer.
- NanoCaptureLigands can be regenerated at least 10 times with Regeneration buffer with minimal loss of binding efficiency.
- Nano-CaptureLigands carry a His-tag; thus, avoid the use of anti-His primary antibodies.
- Briefly centrifuge the Nano-CaptureLigand solution before use.

Protocol

1. Baseline 1:

- Incubate the biosensors for 60 s in 1x Kinetics buffer.

2. Loading

- Dilute the Nano-CaptureLigand to a concentration of 1 µg/mL in 200 µL 1x Kinetics buffer.
- Load the diluted Nano-CaptureLigand onto the biosensors for 60-120 s until a loading response of 1 nm is reached.

Optional: Use the threshold limit function in the FortéBio Data Acquisition software.

3. Quenching (*optional*)

- Incubate the biosensors for 60 s with biocytin (10 µg/mL in 1x Kinetics buffer) .

4. Baseline 2

- Incubate the biosensors for 120 s in 1x Kinetics buffer.

5. Activation

- Activate the biosensors for 120-180 s with the antibody (20 nM in 1x Kinetics buffer).

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

6. Baseline 3

- Incubate the biosensors for 120 s in 1x Kinetics buffer.

7. Association

- Bind different antigen concentrations in 1x Kinetics buffer for 120-600 s.

Note: As a start, use 0.1-250 µg/mL or 1/10-10x K_d of antigen.

8. Dissociation

- Incubate the biosensors for 60-800 s in 1x Kinetics buffer.


Note: Use the wells from step 6, Baseline 3.

Note: Duration of dissociation step depends on the affinity of the analyzed interaction.

9. Regeneration (optional)

- Regenerate the biosensors for 5 s with Regeneration buffer.
- Incubate in 1x Kinetics buffer for neutralization.
- Repeat regeneration 2 times.

Application examples

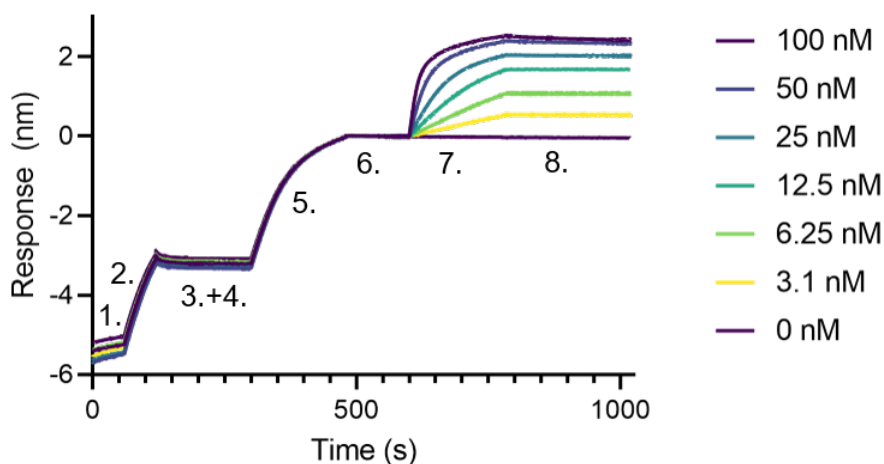


BLI binding kinetics of a human IgG1 anti-IgE antibody to human IgE. A human IgG1 anti-IgE antibody was immobilized using Nano-CaptureLigand human IgG/rabbit IgG, Fc-specific VHH, biotinylated on FortéBio Streptavidin (SA) Biosensors and assayed with different concentrations of human IgE.


Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

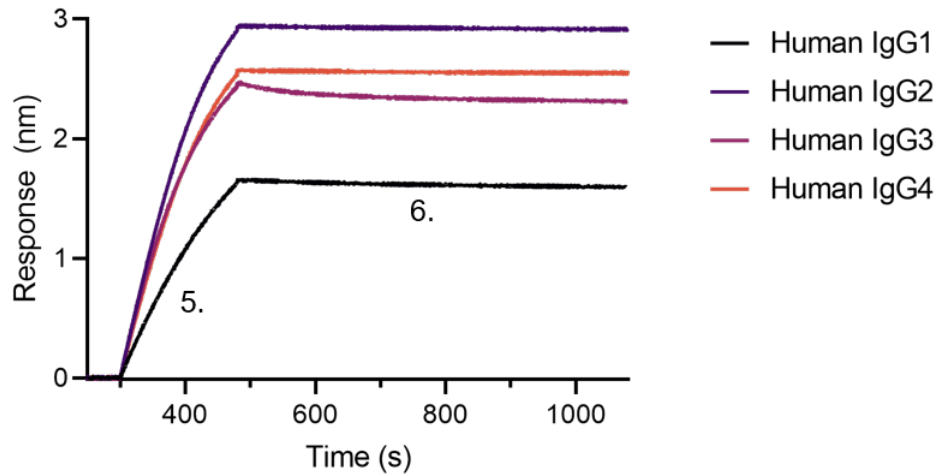

BLI binding kinetics of different human IgG antibodies. Different monoclonal human IgG antibodies were immobilized using Nano-CaptureLigand human IgG/rabbit IgG, Fc-specific VHH, biotinylated on FortéBio Streptavidin (SA) Biosensors. All four human IgG subclasses are stably captured by the Nano-CaptureLigand with negligible dissociation.

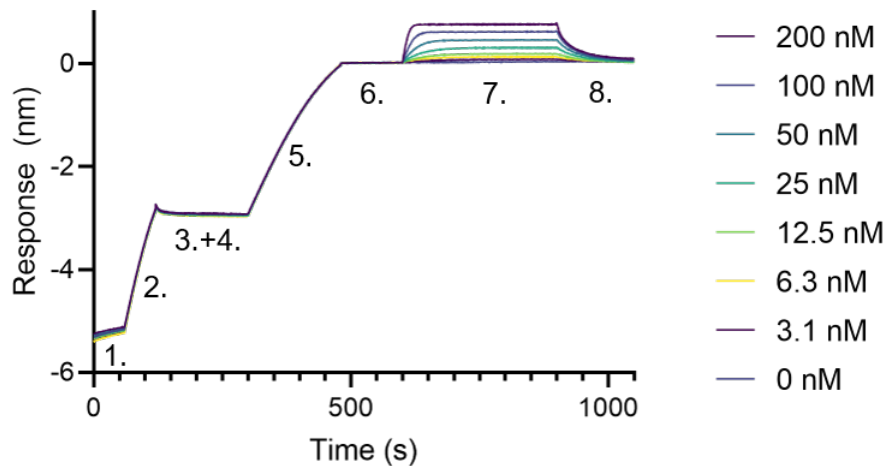


BLI binding kinetics of a monoclonal rabbit IgG anti-DYKDDDDK-tag antibody to a DYKDDDDK-tagged protein. A monoclonal rabbit IgG anti-DYKDDDDK-tag antibody was immobilized using Nano-CaptureLigand human IgG/rabbit IgG, Fc-specific VHH, biotinylated on FortéBio Streptavidin (SA) Biosensors and assayed with different concentrations of a DYKDDDDK-tagged protein.

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

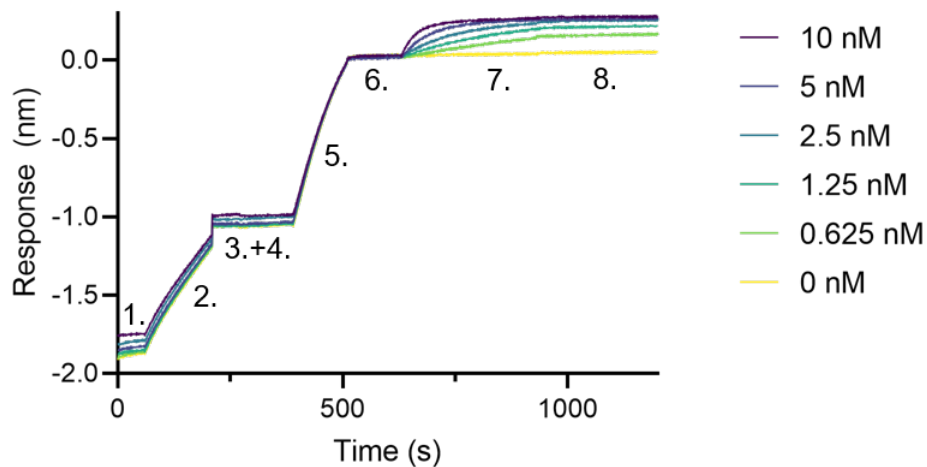

BLI binding kinetics of a polyclonal rabbit IgG anti-GFP antibody to eGFP. A polyclonal rabbit IgG anti-GFP antibody (ChromoTek GFP antibody [PABG1]) was immobilized using Nano-CaptureLigand human IgG/rabbit IgG, Fc-specific VHH, biotinylated on FortéBio Streptavidin (SA) Biosensors and assayed with different concentrations of eGFP (ChromoTek EGFP, recombinant protein).

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

Sandwich ELISA protocol

General notes

- The following protocol was designed for a standard sandwich ELISA. Other types of ELISA or your specific research question may require optimization of particular parameters.
- Use the recommended materials or their equivalents.
- In this protocol, MaxiSorp plates (e.g. Thermo Scientific™ White and Black 384-Well Immuno Plates, #460518) are used that must be coated with avidin or streptavidin first. Alternatively, pre-coated avidin/streptavidin plates can be used.
- Nano-CaptureLigands carry a His-tag; thus, avoid the use of anti-His primary antibodies.
- Recommended volumes for 96-well and 384-well microplates:

Protocol steps	96-well microplate	384 -well microplates
Coating, antigen binding, antibody binding	100 µL	20 µL
Washing, blocking	300 µL	90 µL

- Briefly centrifuge the Nano-CaptureLigand solution before use.

Protocol

1. Avidin coating (*optional*)

- Coat each well of a MaxiSorp plate with 10 µg/mL avidin in PBS at +4°C overnight.
- Wash each well twice with PBS.

2. Blocking

- Block each well with Blocking buffer for 1-2 h at room temperature.
- Wash each well 3 times with Wash buffer.

3. Nano-CaptureLigand coating

- Add 50 nM Nano-CaptureLigand (diluted in Dilution buffer) to each well.
- Incubate for 1 h at room temperature.
- Wash each well 5 times with Wash buffer.

4. Immobilization of capture antibody

- Add the capture antibody (diluted in Dilution buffer) to each well.
- Incubate for 1 h at room temperature.
- Wash each well 5 times with Wash buffer.

Note: Test different concentrations of the capture antibody in an initial experiment.

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

5. Antigen binding

- Add the antigen to each well.
- Incubate for 1 h at room temperature.
- Wash each well 5 times with Wash buffer.

Note: Test different concentrations of the antigen.

6. Binding of primary antibody

- Add the primary antibody to each well and incubate.
- Wash each well 5 times with Wash buffer.

Note: Dilute and incubate the primary antibody as indicated in the manufacturer's manual.

7. Binding of secondary / detection antibody

- Add the secondary / detection antibody to each well and incubate.
- Wash each well 5 times with Wash buffer.

Note: Dilute and incubate the secondary / detection antibody as indicated in the manufacturer's manual.

8. Detection


- Add the appropriate ELISA substrate solution to each well and incubate as indicated in the manufacturer's manual.
- Analyze with a microplate reader.

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

Application examples

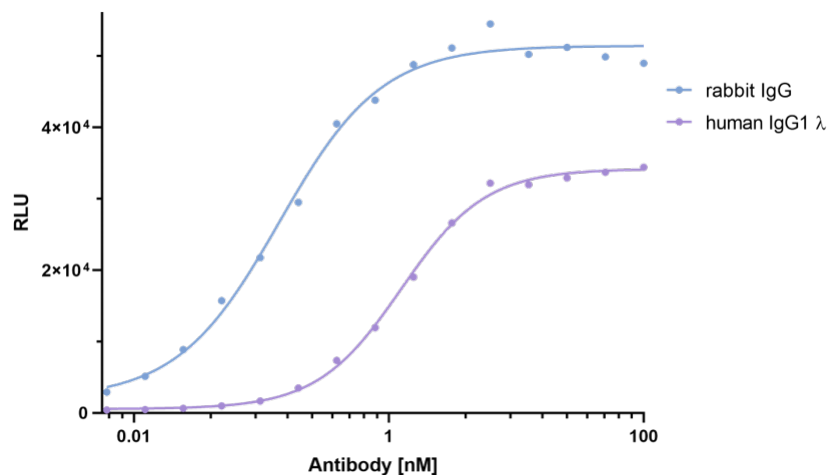

ELISA capture of a human IgG1 lambda and a rabbit IgG antibody using Nano-CaptureLigand human IgG/rabbit IgG, Fc-specific VHH, biotinylated. 50 nM Nano-CaptureLigand human IgG/rabbit IgG, Fc-specific VHH, biotinylated was used for coating on an avidin-coated MaxiSorp plate. Human IgG1 lambda antibody and rabbit IgG antibody were titrated in a 1:2 dilution series and detected with an alkaline phosphatase-conjugated detection antibody.

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

Product overview and related products

Product	Product code
Nano-CaptureLigand™ human IgG/rabbit IgG, Fc-specific VHH, biotinylated	shurbGB-1-10; -100
Nano-CaptureLigand™ human Ig, lambda-LC-specific VHH, biotinylated	shuLB-1-10; -100
Nano-CaptureLigand™ human IgE, VHH, biotinylated	shuEB-1-10; -100
Nano-CaptureLigand™ mouse IgG1, Fc-specific VHH, biotinylated	smsG1B-1-10; -100
Nano-CaptureLigand™ mouse IgG2a, Fc-specific VHH, biotinylated	smsG2aB-1-10; -100
Nano-CaptureLigand™ mouse IgG2b, Fc-specific VHH, biotinylated	smsG2bB-1-10; -100
Nano-CaptureLigand™ mouse IgE, VHH, biotinylated	smsEB-1-10; -100
GFP VHH, biotinylated recombinant binding protein	gtb-250

For product details, information, and ordering visit www.chromotek.com.

Nano-CaptureLigand™ human IgG/ rabbit IgG, Fc-specific VHH, biotinylated



Product code: shurbGB-1

Contact

support@chromotek.com

ChromoTek GmbH
Am Klopferspitz 19
82152 Planegg-Martinsried
Germany
phone: +49 89 124 148 80
fax: +49 89 124 148 811

ChromoTek Inc.
62-64 Enter Lane
Islandia, NY 11749
USA
phone: 631 501 1058
fax: 631 501 1060

Disclaimer

Only for research applications, not for diagnostic or therapeutic use!

ChromoTek and GFP-Trap®, RFP-Trap®, Myc-Trap®, Spot-Trap®, Spot-Tag®, Spot-Label®, Spot-Cap®, Nano-Secondary®, F2H® Kit, and Chromobody® are registered trademarks and Nano-CaptureLigand™ is a trademark of ChromoTek GmbH. Nanobody is a registered trademark of Ablynx, a Sanofi company. Octet® is a registered trademark of ForéBio, a Sartorius brand. Dynabeads™ is a trademark of Life Technologies AS, a part of Thermo Fisher Scientific Inc. Alexa Fluor® is a registered trademark of Life Technologies Corporation, a part of Thermo Fisher Scientific Inc. Other suppliers' products may be trademarks or registered trademarks of the corresponding supplier each. Statements on other suppliers' products are given according to our best knowledge.