

Buffer and Solution Preparation

Instructions

1. Prepare fresh Transcription Factor Fixation/Permeabilization working solution by diluting Transcription Factor Fixation/Permeabilization Concentrate (1 part) with Transcription Factor Fixation/Permeabilization Diluent (3 parts). You will need 1 mL of the Transcription Factor Fixation/Permeabilization working solution for each sample.
2. Prepare a 1X working solution of Flow Cytometry Perm Buffer by diluting the 10X concentrate with distilled water prior to use. You will need 3-5 mL of Flow Cytometry Perm Buffer working solution for each sample.

More guides available upon request at ptglab.com