

SignalBright Plus Chemiluminescent Substrate

Contents:

- 1x Chemiluminescent Substrate Solution, 1x Stable Peroxide Solution
- Recommended primary antibody concentration: **0.02-0.5 µg/ml**
- Recommended HRP-conjugated secondary antibody concentration: **2-20 ng/ml**

Instructions

- Add equal parts of the Chemiluminescent Substrate Solution and the Stable Peroxide Solution to create a working solution
- Create enough working solution with to allow a minimum of 0.1ml per membrane cm².
- Allow the working solution to come to room temperature in complete darkness
- Apply the working solution to the membrane and allow to develop for 5 minutes

More guides available upon request at ptglab.com