

# PURIFICATION OF PROTEINS FROM INCLUSION BODIES

[www.ptglab.com](http://www.ptglab.com)

Recipes for all solutions (highlighted) in **bold** are included at the end of the protocol.

1.
  - a. Suspend the cell pellet (from 1 L culture) in 30–35 ml of **PBST buffer**.
  - b. Sonicate cells in an ice-bath at 200 W for 8 min.
  - c. Centrifuge cell lysate for approximately 15 min at 8000 rpm, 4°C. Discard the supernatant.
  - d. Re-suspend the pellets in 30 ml **TNMFx-2M Urea buffer** before transferring to a 50 ml centrifuge tube. Rotate for 30 min at 4°C.
  - e. Sonicate the solution in an ice-bath at 200 W for 1–4 min.
  - f. Rotate for 30 min at 4°C.
  - g. Centrifuge for 15 min at 8000 rpm, 4°C. Discard the supernatant.
  - h. Re-suspend the pellets in 30ml of **TNMFx-0.1% Triton-X100**. Rotate for 30 min. at 4°C.
  - i. Centrifuge for 15 min at 8000 rpm, 4°C. Discard the supernatant. Vortex and wash the pellets with 2x volumes of dH<sub>2</sub>O before transferring to a 10 ml centrifuge tube. Centrifuge at 8000 rpm for 10 min.
  - j. Repeat washing until the supernatant becomes clear. Collect the pellets.
  - k. Repeat washing until the supernatant becomes clear. Collect the pellets.
  - l. Dissolve the proteins depending on intended application:
    - For immunization, dissolve in 1.5x volumes of **8 M urea** (pH 8.0)
    - For antibody purification, incubate in 2x volumes of **PBS with 2% Sarkosyl** overnight at 4°C. Collect the supernatant by centrifugation at 12000 rpm for 15 min.

## Solutions

TNMFx-2M Urea	For 1000 ml
50 mM Tris-base	6.06 g
150 mM NaCl	8.77 g
1 mM EDTA	0.37 g
2 M Urea	120.20 g
Adjust to pH 8.0	
Add ddH <sub>2</sub> O to 1000 ml	

## Contact

**US Head Office**  
proteintech@ptglab.com

**United Kingdom**  
europe@ptglab.com

**Proteintech China**  
proteintech-cn@ptglab.com

# PURIFICATION OF PROTEINS FROM INCLUSION BODIES

[www.ptglab.com](http://www.ptglab.com)

## Solutions

PBST buffer	For 1000 ml
58 mM Na <sub>2</sub> HPO <sub>4</sub> • 12H <sub>2</sub> O	20.77 g
17 mM NaH <sub>2</sub> PO <sub>4</sub> • 2H <sub>2</sub> O	2.65 g
68 mM NaCl	3.98 g
1% Triton-X100	10 ml
Adjust to pH 7.4	
Add ddH <sub>2</sub> O to 1000 ml	

TNMFx-0.1% Triton X100	For 1000 ml
50 mM Tris	6.06 g
150 mM NaCl	8.8 g
1 mM EDTA	0.4 g
0.1% Triton-X100	1 ml
Adjust to pH 8.0	
Add ddH <sub>2</sub> O to 1000 ml	

PBS with 2% Sarkosyl	For 200 ml
58 mM Na <sub>2</sub> HPO <sub>4</sub> • 12H <sub>2</sub> O	4.19 g
17 mM NaH <sub>2</sub> PO <sub>4</sub> • 2H <sub>2</sub> O	0.53 g
68 mM NaCl	0.80 g
2% Sarkosyl	4.00 g
Adjust to pH 8.0	
Add ddH <sub>2</sub> O to 200 ml	

8 M Urea	For 200 ml
Urea	96.08 g
Adjust to pH 8.0	
Add ddH <sub>2</sub> O to 200 ml	